



BROWN
Division of Biology
and Medicine

REES Monitoring System

The REES monitoring system can be accessed by both the REES web log-in, and by Phone

The REES web log-in is: <https://preescit1.ad.brown.edu/>

The phone numbers for the respective nodes are as following:

- Node 1 – 401-863-1651
- Node 2 – 401-863-9135
- Node 3 – 401-863-6523

To access the REES web log in:

- Open up the provided link in your web browser
- Type in your user name and password (If you do not have a user name or password, or if you do not remember it, please contact the office of Bio Med FPO at ext. 3-2483)
- To see your respective inputs, click on the select a node tab, and select the node for your respective equipment. (If you do not know what node your inputs are on, please contact the office of BioMed FPO at ext. 3-2483)
- Once you see the input numbers on the screen, bringing the mouse over the input number will allow you to see the current state of the input in the box at the bottom of the screen.
- Clicking on the respective input will allow you to do the following: Inhibit, or deactivate the alarm, or reset the input. Resetting the input will typically be used to cancel out an alarm inhibit.
- To see a graph of the input data, click on the blue tab “Graph Input Readings” on the left of the screen. You will be able to select the input you would like to graph, as well as determine the window

To access the REES system by Phone:

- Call the number for you respective node (If you do not know what node your inputs are on, please contact the office of Bio Med FPO at ext. 3-2483)
- The system will answer and require you to enter your user 4 digit ID followed by the # (If you do not know what node your inputs are on, please contact the office of Bio Med FPO at ext. 3-2483)



BROWN

Division of Biology and Medicine

- The phone will give you 5 options:
- “Press 1 to check global status” –
- “Press 2 to check single input status” - This will allow you to get the current reading of the input
- “Press 3 to inhibit an input” – This will allow you to inhibit, or silence the alarm, for the input for a duration up to 9999 hours
- “Press 4 to enable an input” – This will allow you to reset the input if it is currently inhibited
- “Press 5 to hang up”

The Rees Monitoring system will contact the lab by both email and by phone if an input is in alarm. When an input is in alarm for 15 minutes, the system will send out an email to the people on the email list determined by the lab. The system has the capability for up to three people to be on the email list.

When an input is in alarm for 25 minutes, the system will begin to call the phone list provided to our office by each respective lab. When the system calls, the five options will be the exact same as if you were to call the system. The only difference is that the system will tell you what input is in alarm, prior to allowing you to enter your telephone ID.

If you have any questions on your respective labs existing email or phone list, or if you would like to make any changes to the existing lists, please send your updated list to our office at biomed_facilities@brown.edu.