

Transferring files from nmr computers to christie:

1. Double click on SSH Secure File Transfer icon.
2. Click quick connect.
3. Type in nmr computer <name>.chem.brown.edu where <name> is surya, oksana, katarina, or tara. Enter your login (user name). Click connect. Enter password.
4. Find your files on the nmr computer on the left side of the screen under <login> nmr where <login> is your nmr login. Click nmr. Click on the files you want to transfer on the right side of the screen.
5. Open another window by double clicking My Computer icon.
6. Double click local disk C. Double click Bruker, double click Topspin, double click data, double click guest, double click nmr. This is the Topspin directory where files should be transferred. Drag the files from the first window to the second window.
7. Close the second window under file. Disconnect the first window under file. Click yes. Exit under file.